



# **INSTRUCTION MANUAL FOR CHIRALPAK® MA(+)**

# Please read this instruction sheet completely before using this column

**Column Description** 

Packing composition:

N,N-dioctyl-L-alanine coated on **3µm silica-gel**.



Column size:

50mm L x 4.6mm i.d.

Shipping solvent:

The column is filled with  $2mM CuSO_4$  aq. solution.

All columns have been pre-tested before packaging. Test parameters and results, as well as the Column Lot Number, are included on a separate (enclosed) page.

### CAUTION

This column contains an amino acid derivative coated on silica-gel as the packing material. The entire HPLC system including the injector and the injection loop must be flushed with a mobile phase compatible with the column and its storage solvent prior to connecting.

Do not use mobile phases other than those listed in this document. Otherwise, the column may be damaged by dissolving the amino acid derivative. Strong-basic compounds are likely to deteriorate the silica-gel used in this column.

# **Standard Analytical Conditions**

Mobile phase	: $2mM$ CuSO <sub>4</sub> aq. solution
Operating temperature	: 0 - 40°C
Column pressure	: Max. 150kg/cm <sup>2</sup>
Flow rate	: 1.0mL/min
Flow direction	: As indicated on column label

# **Mobile Phases**

Suitable mobile phases :	
--------------------------	--

 $CuSO_4$  aq. solution  $CuSO_4$  aq. solution / methanol  $CuSO_4$  aq. solution / acetonitrile = 0.1 - 2.0mM = 100/0 - 85/15 (v/v) = 100/0 - 85/15 (v/v)

#### NOTES:

- $\Rightarrow$  Retention time is generally shorter with acetonitrile than with methanol and with a higher content of organic solvent.
- $\Rightarrow$  Eluent solution should be filtered through a 0.45 mm membrane filter.
- $\Rightarrow$  High content of organic solvent may cause precipitation of CuSO<sub>4</sub> which makes the tubing clogged.

#### Sample

The sample solution should be filtered through a 0.45 mm membrane filter as a rule.

#### **Column Cleaning**

If the column has been contaminated, wash the column with 15% acetonitrile in  $H_2O(v/v)$  at a flow rate of 0.5mL/min for 3 hours.

#### Storage

The column can be stored at room temperature. When the column is not in use, it is recommended to fill it with a 2mM  $CuSO_4$  aq. solution.

#### **Important Notice**

 $\Rightarrow$  STRONGLY BASIC solvent modifiers or sample solutions MUST BE AVOIDED, because they are likely to damage the silica gel used in this column.

#### Operating this column in accordance with the guidelines outlined here will result in a long column life.

⇒ If you have any questions about the use of these columns, or encounter a problem, contact:

In the USA: <u>questions@chiraltech.com</u> or call 800-6-CHIRAL In the EU: <u>cte@chiral.fr</u> or call +33 (0)3 88 79 52 00 In India: <u>chiral@chiral.daicel.com</u> or call +91-40-2338-3700

#### Locations:

#### North/Latin America

Chiral Technologies. Inc. 800 North Five Points Road West Chester, PA 19380 800 6 CHIRAL Tel: 610-594-2100 Fax: 610-594-2325 chiral@chiraltech.com www.chiraltech.com **Europe** Chiral Technologies Europe Parc d'Innovation Bd Gonthier d'Andernach 67400 Illkirch Cedex, France Tel: +33-388-795-200 Fax: +33-388-667-166 cte@chiral.fr www.chiral.fr India

Daicel Chiral Technologies (India) Pvt. Ltd. Lab No. 4A, Phase III IKP Knowledge Park Genome Valley, Turkapally, Shameerpet, Ranga Reddy Dist. Hyderabad-500 078, Telangana Tel: +91-40-2338-3700 Fax: +91-40-2348-0104 chiral@chiral.daicel.com

CHIRALCEL, CHIRALPAK and CROWNPAK are registered trademarks of DAICEL CORPORATION